Check Stab α 2016 iLife

Caratteristiche tecniche - Technical characteristics

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Data given in this brochure are indicative and can vary at the discretion of the manufacturer at any time and without obligation or notice.

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Functions of the Check Stab α2016 iLife

- *Quick and easy measurement of tartaric acid stability and calcium saturation using electrochemical conductivity measurements at various conditions.*
- *Mini Contact: Determines KHT precipitation expressed in μSiemens and %. This value indicates the relative stability of the wine sample being tested.*
- *Mini contact with forecast: Determines KHT precipitation expressed in μSiemens and %. It uses CheckFuture® to predict future knock down of crystals.*
- *Saturation Point: Automatically, the analysis calculates the saturation temperature of the test sample expressed in °C. It also has the possibility to evaluate the data with the Total Polyphenols Index.*
- *Saturation Point of Calcium: This analysis calculates automatically the value of saturation temperature of calcium.*
- *Special Functions: ISS®, ISTC50.*

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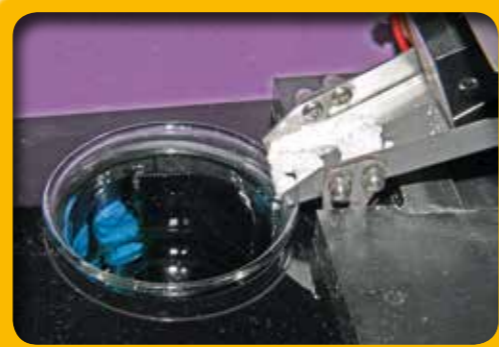
Some characteristics of the Check Stab α2016 iLife

The operation and management of data is done with a PC with a Windows operating system. The instruments can be connected to the PC by LAN. Possibility to connect to Bar Code reader to insert sample data. Possibility to connect more than one instrument together (max 16 units), with each unit doing an independent analysis. Automatic dispensing system for adding KHT without the intervention of the lab technician. Stand by function for energy savings.

Due to numerous requests from many of our customers, **CheckStab Instruments**®, leader by virtue of new technologies and experimentation, has developed a new innovative instrument for the analysis of tartaric stability in wines. The instrument is equipped with a 9,9" touch screen colour display, interfaced with CheckStab.NET® software by means of LAN, which means it can also be connected directly to the winery's computer network. **Check Stab α2016 iLife** is a completely automatic analyzer capable of satisfying all types of wineries, small or large. The laboratory can connect from 1 to 16 instruments together, all controlled by one software CheckStab Net®, with the data stored on one single Data Base. The instrument has an automatic KHT dispensing system.

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View of the automatic KHT dispenser

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Acquiring and recording measurement data is becoming more and more important. **CheckStab Instruments**® has developed a new analytical software system to manage data. This permits the lab to study, store, and comment on an analysis with extreme simplicity. The entire system is managed by use of a personal computer.

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What Check Stab α2016 iLife and Check Stab.NET can do:

- *Contemporaneous management of connected instruments*
- *Possibility to do at the same time several analysis, with independent parameter research*
- *Visualize measurement data [conductivity, wine temperature, bath temperature, pre-set temperature, etc.]*
- *Visualize graphics in real time*
- *Possibility to start the analysis directly from instrument and automatic synchronization with the PC.*
- *Visualize user status, using incorporated graphic symbols (State of cooler, heater, probe positions, etc.).*
- *Memorize the analysis on the data-base, with the possibility to export to a text file.*
- *Possibility to get from Check Stab analyses done, to visualize results in graphic form, to visualize results in text form, to save on a disk, export reports in various formats (PDF, Excel, Text) or print the graphics using a printer.*
- *Visualize the saturation analysis in a normal graphics mode or in a differential mode.*
- *Visualize the mini contact analysis in absolute mode or in relative mode.*
- *Visualize the analysis in single mode or multiple mode, up to a maximum of 10 analysis contemporaneously.*
- *Interface for LIMS, sends reports automatically by e-mail.*